

### Compact NIR MCP-PMT Series Featuring with Fast Time Response

#### FEATURES

- High Speed  
Rise Time: 170 ps (Typ.)  
T.T.S. (Transit Time Spread):  $\leq 70$  ps (FWHM)<sup>Ⓐ</sup> (Typ.)

#### APPLICATIONS

- Fluorescence, Fluorescence Lifetime
- Photoluminescence
- LIDAR



▲ H10840-68/-69



▲ NIR MCP-PMT R3809U-68/-69  
Included in H10840-68-69

Figure 1: Spectral Response


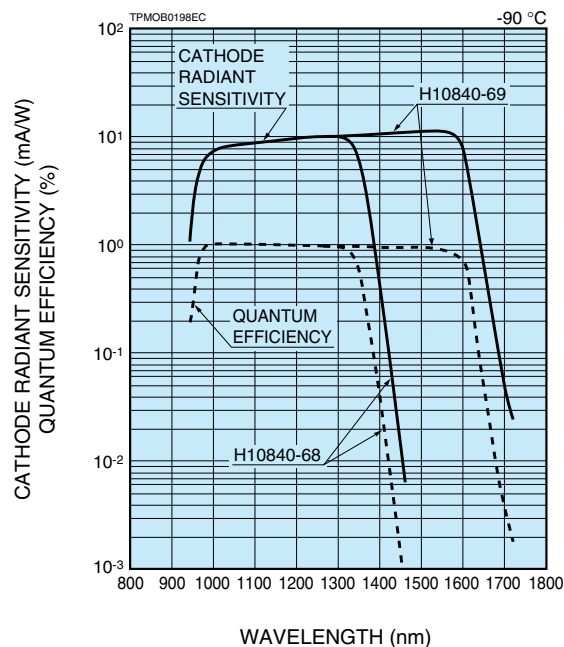


Figure 2: Typical Gain

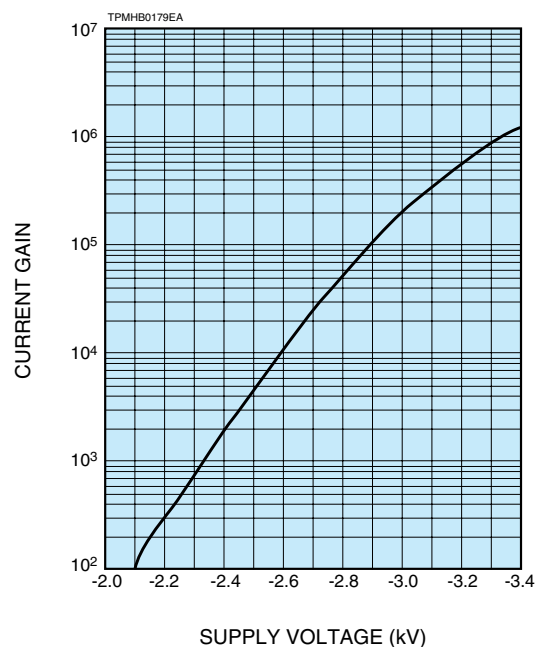

Subject to local technical requirements and regulations, availability of products included in this promotional material may vary. Please consult with our sales office. Information furnished by HAMAMATSU is believed to be reliable. However, no responsibility is assumed for possible inaccuracies or omissions. Specifications are subject to change without notice. No patent rights are granted to any of the circuits described herein. ©2009 Hamamatsu Photonics K.K.

# NIR MCP-PMT WITH COOLER H10840-68/-69

## SPECIFICATIONS

### GENERAL

Parameter	H10840-68	H10840-69	Unit
Spectral Response	950 to 1400	950 to 1700	nm
Photocathode Material	InP / InGaAsP	InP / InGaAs	—
Window Material	Borosilicate Glass		—
Effective Area of PMT	φ2		mm
Stage of MCP <sup>®</sup>	2		—
Operating Ambient Temperature	0 to 40		°C
Storage Temperature	-20 to +50		°C

### MAXIMUM RATING

Parameter	H10840-68	H10840-69	Unit
PMT Supply Voltage	-3400		V
Average PMT Anode Current	20		nA

### CHARACTERISTICS (at -3000 V, -90 °C)

Parameter		H10840-68			H10840-69			Unit
		Min.	Typ.	Max.	Min.	Typ.	Max.	
Cathode Sensitivity <sup>©</sup>	Quantum Efficiency	0.1	1.0	—	0.1	1.0	—	%
	Radiant	1.0	10.5	—	1.2	12.1	—	mA/W
Gain		1 × 10 <sup>5</sup>	2 × 10 <sup>5</sup>	—	1 × 10 <sup>5</sup>	2 × 10 <sup>5</sup>	—	—
Anode Dark Count <sup>®</sup>		—	2 × 10 <sup>4</sup>	1 × 10 <sup>5</sup>	—	5 × 10 <sup>4</sup>	2.5 × 10 <sup>5</sup>	s <sup>-1</sup>
Voltage Driver Current		—	—	100	—	—	100	μA
Time Response	Rise Time <sup>®</sup>	—	170	—	—	170	—	ps
	Fall Time <sup>®</sup>	—	450	—	—	450	—	ps
	Transit Time Spread <sup>®</sup>	—	70	100	—	70	100	ps

<sup>®</sup> Transit-time spread (T.T.S.) is the fluctuation in transit time between individual pulse and specified as an FWHM (full width at half maximum) with the incident light having a single photoelectron state.

<sup>®</sup> Two microchannel plates (MCP) are incorporated as a standard but we can provide it with either one or three MCPs as an option depending upon your request.

<sup>©</sup> At 1300 nm (H10840-68), at 1500 nm (H10840-69)

<sup>®</sup> At 30 minutes after high voltage is applied with shutter closed

<sup>®</sup> This is the mean time difference between the 10 % and 90 % amplitude points on the output waveform for full cathode illumination.

<sup>®</sup> This is the mean time difference between the 90 % and 10 % amplitude points on the tailing edge of the output waveform for full cathode illumination.

<sup>®</sup> I.R.F. stands for Instrument Response Function which is a convolution of the δ pulse function (H(t)) of the measuring system and the excitation function (E(t)) of a laser. The I.R.F. is given by the following formula:

$$I.R.F. = H(t) \times E(t)$$

We specify the I.R.F. as an FWHM of the time distribution taken by using the measuring system in Figure 6 that is Hamamatsu standard I.R.F. measurement. It can be temporary estimated by the following equation:

$$(I.R.F. (FWHM))^2 = (T.T.S.)^2 + (T_w)^2 + (T_j)^2$$

where  $T_w$  is the pulse width of the laser used and  $T_j$  is the time jitter of all equipments used. An I.R.F. data is provided with the tube purchased as a standard.

### COOLER

Parameter	Description / Value	Unit
Coolant Medium	Liquid Nitrogen (LN <sub>2</sub> )	—
Temperature Controllable Range	0 to -100 (continuously adjustable)	°C
Cool-down Time	About 2 (-90 °C setting)	h
Liquid Nitrogen Consumption	Approx. 0.75 (-100 °C setting)	L/h
Dry Nitrogen	Gas Pressure	35
	Consumption	47 L (14.7 MPa) / 10 h
Holder	-HV Connector	SHV-R
	Signal Connector	SMA-R
	Load Resistor	Open
Power Consumption	15	VA
Operating Ambient Temperature	Less than +40	°C
Weight	Cooling Unit	Approx. 6
	Controller, etc.	Approx. 11
System Configuration	Cooling Unit, Controller, Control Cable, Solenoid Control Cable, Rubber Tube, Insulated Transfer Hose, LN <sub>2</sub> Transfer Head	—

Figure 3: Typical Output Waveform


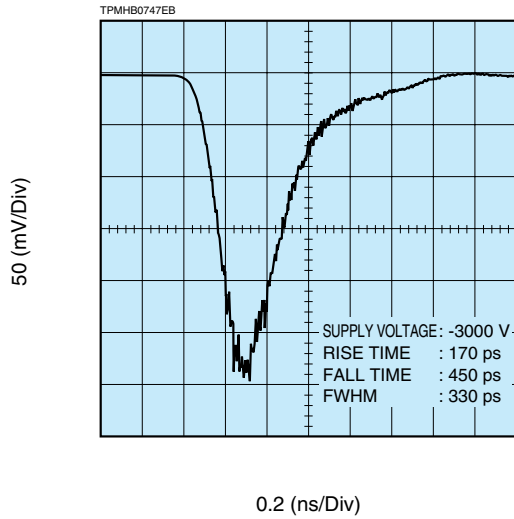


Figure 4: Block Diagram of Output Waveform Measuring System

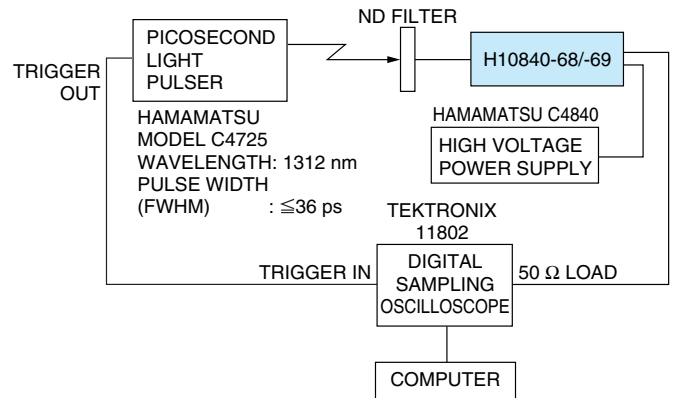

TPMHC0230ED

Figure 5: Typical Instrument Response Function (I.R.F.)<sup>®</sup>


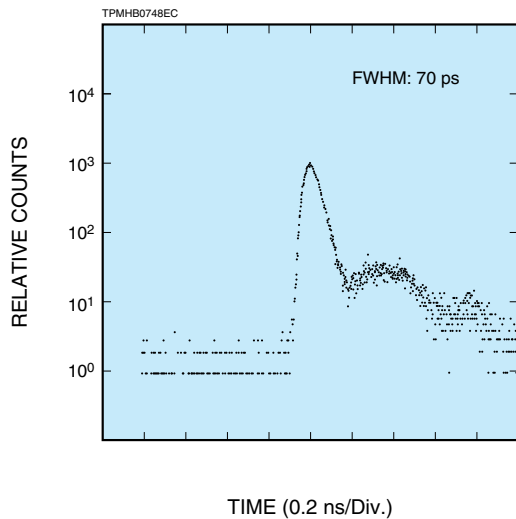


Figure 6: Block Diagram of I.R.F. Measuring System

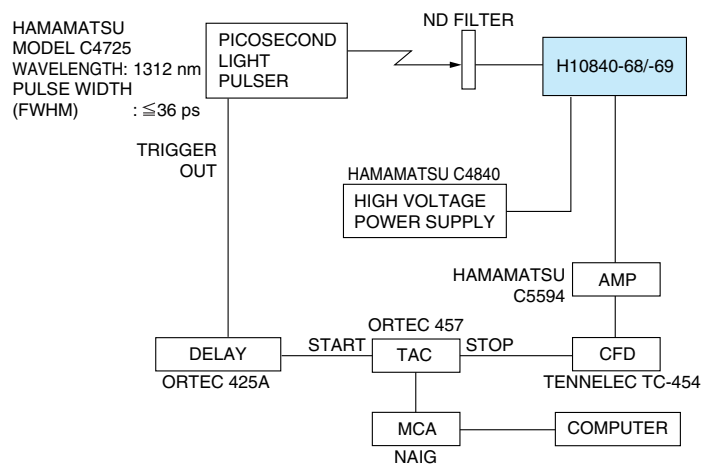


TPMHC0231ED

# NIR MCP-PMT WITH COOLER H10840-68/-69

Figure 7: Dimensional Outlines (Unit: mm)

## ●COOLING UNIT

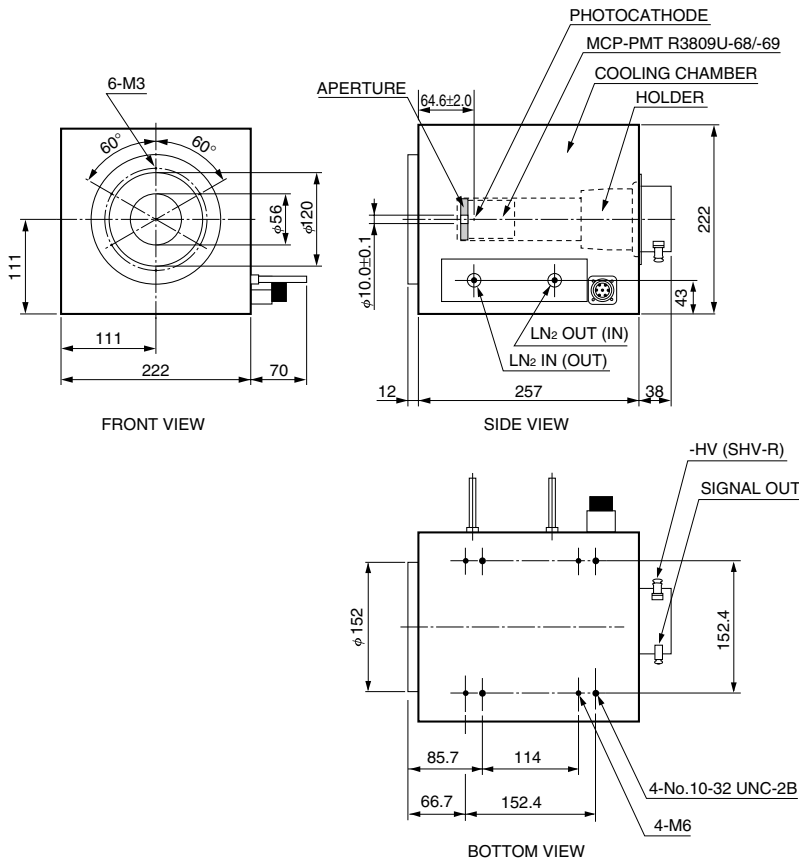

## ●CONTROLLER

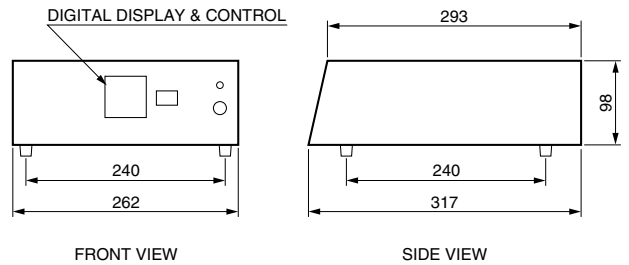


Figure 8: System Configuration

### \*Other accessories required (Not supplied)

#### ●LN<sub>2</sub> dewar


Non-pressurized dewar having a capacity of 10 to 50 liters, and the neck outer diameter between 35 and 40 mm.

#### ●High voltage power supply for the photomultiplier tube (negative high voltage)

Output voltage: more than -3400 V  
Output current: more than 0.2 mA  
Low ripple, High stability

#### ●Dry nitrogen gas, pressure regulator (secondary pressure 35 kPa)

In order to supply a proper amount of liquid nitrogen to the cooling unit, an external pressure needs to be added to the dewar. A pressure regulator capable of reducing a secondary pressure to 35 kPa is necessary when used with a dry nitrogen gas container. Connect the 3/8" rubber tube to the exit of the pressure regulator.

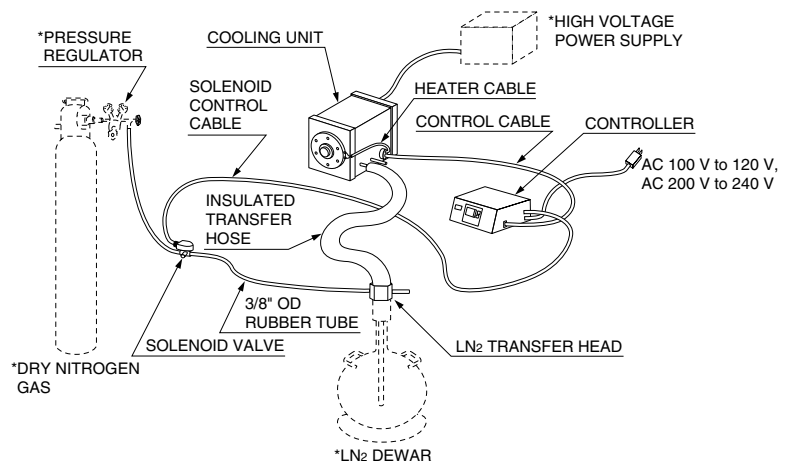

# HAMAMATSU

WEB SITE [www.hamamatsu.com](http://www.hamamatsu.com)

HAMAMATSU PHOTONICS K.K., Electron Tube Division

314-5, Shimokanzo, Iwata City, Shizuoka Pref., 438-0193, Japan, Telephone: (81)539/62-5248, Fax: (81)539/62-2205

U.S.A.: Hamamatsu Corporation: 360 Foothill Road, P. O. Box 6910, Bridgewater, N.J. 08807-0910, U.S.A., Telephone: (1)908-231-0960, Fax: (1)908-231-1218 E-mail: [usa@hamamatsu.com](mailto:usa@hamamatsu.com)

Germany: Hamamatsu Photonics Deutschland GmbH: Arzbergerstr. 10, D-82211 Herrsching am Ammersee, Germany, Telephone: (49)8152-375-0, Fax: (49)8152-2658 E-mail: [info@hamamatsu.de](mailto:info@hamamatsu.de)

France: Hamamatsu Photonics France S.A.R.L.: 19, Rue du Saule Trapu, Parc du Moulin de Massy, 91882 Massy Cedex, France, Telephone: (33)1 69 53 71 00, Fax: (33)1 69 53 71 10 E-mail: [infos@hamamatsu.fr](mailto:infos@hamamatsu.fr)

United Kingdom: Hamamatsu Photonics UK Limited: 2 Howard Court, 10 Tewin Road Welwyn Garden City Hertfordshire AL7 1BW, United Kingdom, Telephone: 44-(0)1707-294888, Fax: 44(0)1707-325777 E-mail: [info@hamamatsu.co.uk](mailto:info@hamamatsu.co.uk)

North Europe: Hamamatsu Photonics Norden AB: Smidesvägen 12, SE-171-41 SOLNA, Sweden, Telephone: (46)8-509-031-00, Fax: (46)8-509-031-01 E-mail: [info@hamamatsu.se](mailto:info@hamamatsu.se)

Italy: Hamamatsu Photonics Italia: S.R.L.: Strada della Moia, 1/E, 20020 Arese, (Milano), Italy, Telephone: (39)02-935 81 733, Fax: (39)02-935 81 741 E-mail: [info@hamamatsu.it](mailto:info@hamamatsu.it)

TACCC0090EE

TPMH1312E02  
MAY 2009 IP